








ARM – Advanced Real Time Monitor 2011

Icon manual

● Live setting






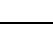
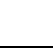
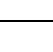

	Start	Start live image
	Stop	Stop live image

● Image storage icon



	Save Document	Save unique ARM images(*.arm).
	Open Document	Open unique ARM images(*.arm). Open general images(Jpg, jpeg, tif, tiff, bmp, png).
	Capture	According to the setting of  (Capture option), save present monitor images.
	Sequency Capture	According to the setting of  (Capture option), save present monitor images with time interval specified.
	Capture option	Set up the folder for Capture images, assign automatic file name, select format of image storage(jpg, tiff, bmp. png), assign storage time interval(minimum time : 2 sec)

ARM – Advanced Real Time Monitor 2011





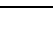



● Measurement icon

	Show only	Click when moving out of measuring state.
	Point	Measure the number of points and coordinates when running ARM at the first.
	Distance	Measure the distance between two points.
	angle	Measure the angle.
	Triangle	Figure out the area of triangle by clicking three points.
	Rect	Figure out the area of square by clicking two points.
	Circle	Calculate the diameter and area of circle by clicking the center point and another one point.
	Circle3d	Calculate the area of circle by clicking three points.
	Poly	Figure out the area by clicking several points.(click with pressing Ctrl, calculation is completed under recognition as the last point.)



● Setting icon

	Calibration	Set the calibration information.(Setting by magnification)
	None	No drawing a base line.

ARM – Advanced Real Time Monitor 2011

	Cross line	Mark cross line in the middle of the image.
	Grid line	Mark grid line at the image.
	Color	Select color for cross line and grid line.
	Sheet	Setting for showing or hiding of excel sheet at the bottom of monitor.
	Ruler position	Select the scale bar location.(None, up-left, bottom-left, up-right, bottom-right)
	Ruler size	Set the scale bar size.
	Zoom	Select the magnification entered in calibration. (Setting by magnification of microscope)
	External tools	Open the outside software.(in case of software using images of camera, check 'Stop Camera Monitoring')

● Excel icon

	Excel	Save as CSV file.
	Clear sheet	Delete all measuring items.

